

## Axio Vert SOP

### Set-up

- To turn on, switch the main power on left side at the back, turn the LED lights on by pressing in the on/off button on the right side at the front and switch from 'eco' to 'permanent' and then adjust the brightness.
- To leave in standby when not in use, switch the bulb from 'permanent' to 'eco'.
- Turn the 'nose' of the microscope to select the objective.
- Turn the turret wheel at the bottom of the microscope to select the reflector module (TL/ GFP/ mRFP).

### Acquisition

- Logon to the computer using your GUID and password.
- Open up the **ZEN blue** software, and select **ZEN Pro** for image acquisition.
- In **live** mode, set the exposure, then select **snap** to capture an image...